

Product Information Sheet

Trade Name(s):	In Vivo Gripper
SKU Number:	MDL-IVG-001
Product description:	Tailored for In Vivo Researchers: Specifically designed for the sensitive nature of in vivo research, ensuring gentle and precise handling of your research subjects. Improves your animal handling skill level, reduces stress on animal, and increases handler safety.

MediLumine Inc.


5795 Av. De Gaspé,
Suite #228
Montreal, QC,
Canada H2S 2X3
Phone: +1 514 929-9744
Fax: +1 514 360-6014
www.medilumine.com

The patented innovative perforated design lets mice instinctively grip the platform, significantly reducing their stress compared to other common methods and making them easier to handle.

Increased Handler Safety: Protect yourself from bites and scratches thanks to a more secure handling platform.

Our platform provides a humane approach to rodent handling, giving you peace of mind to focus on what's most important—your research. Transparent with white opacity and reddish hue under white light illumination due to the scattering of blue light.

Proposed Usage

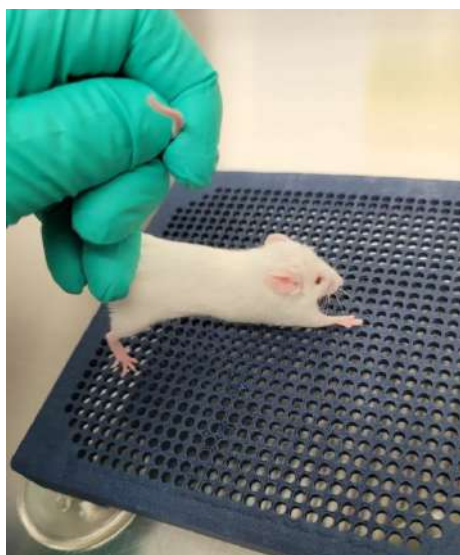

For complete restraint, place the mouse on the in vivo gripper. Securely hold the base of the tail and gently pull back to encourage the animal to grip. Using your other hand, grasp the loose skin at the back of the neck between your thumb and forefinger.